

Index

Sprint ZH 2.1.4_Rev10	2
Improvement	2
Bug	2
New feature	2
Sprint ZH 2.1.4_Rev9	3
New feature	3
Improvement	3
Bug	3
Improvement	3
New feature	3

Sprint ZH 2.1.4_Rev10

Improvement

- [\[ZH-2178\]](#) - Driver TCP Server channel: Limit accepted sockets/connections to 1

Bug

- [\[ZH-2180\]](#) - Driver HL7, HOST_OUT: Sending HL7 results is blocked when an error appears in MeLinDA
- [\[ZH-2183\]](#) - ZH, Daily Work: BlueDiverII - Does not recognize having a pending sample from that machine

New feature

- [\[ZH-2012\]](#) - ZH IFA: Collect and reflect the ZenitPRO warnings regarding some samples

Sprint ZH 2.1.4_Rev9

New feature

- [\[ZH-2170\]](#) - BlueDiverII Driver: Adaptations of the BlueDiver protocol to work with BlueDiverII

Improvement

- [\[ZH-2168\]](#) - Driver HL7, Zenit PRO: Adaptations v1.6.8, change datetime format adding hour minute and second

Bug

- [\[ZH-2147\]](#) - ZH, Configuration, Analyzers, Analyzer / Reagents: A reagent not configured warning is output for an assay not related to the machine
- [\[ZH-2154\]](#) - ZH IFA: Adding a subtest appears without the positivities

Improvement

- [\[ZH-2121\]](#) - ZH, Serumbank, Add tube: Search for the next free position starting from the position occupied by the last added sample
- [\[ZH-2171\]](#) - Driver HL7, ZenitPRO / Fast: Do not process mitosis images
- [\[ZH-2172\]](#) - Driver G-Sight Driver: Receive JP2 images through the WS and not over the network

New feature

- [\[ZH-2025\]](#) - Driver HL7, HOST: Send image analyzer index to LIS