

Index

| | |
|-----------------------|---|
| Sprint ZH 2.1.4 Rev11 | 2 |
| Bug | 2 |
| Improvement | 2 |
| Sprint ZH 2.1.4_Rev10 | 2 |
| Bug | 2 |
| Improvement | 2 |
| New feature | 2 |
| Sprint ZH 2.1.4_Rev9 | 3 |
| Bug | 3 |
| Improvement | 3 |
| New feature | 3 |

Sprint ZH 2.1.4 Rev11

Bug

- [\[ZH-2188\]](#) - ZH IFA: Horizontal Scrollbar insufficient if there are many positivities
- [\[ZH-2190\]](#) - ZH IFA: Certain phantom dilutions are not sent to HOST when end the session
- [\[ZH-2192\]](#) - ZH IFA: PHADIA samples are not displayed in the ZH IFA history
- [\[ZH-2193\]](#) - Driver BlueDiverII: Error when generating session with samples that do not have an associated patient
- [\[ZH-2194\]](#) - Drivers: When installed on a new PC and the IP of the Webservice is configured, the configuration file is saved in the wrong place
- [\[ZH-2195\]](#) - ZH IFA: A warning icon appears when creating phantom dilutions or adding subassays

Improvement

- [\[ZH-2186\]](#) - ZH & ZH IFA & Driver Zenit PRO / Fast: Process and visualize mitosis images

Sprint ZH 2.1.4_Rev10

Bug

- [\[ZH-2180\]](#) - Driver HL7, HOST_OUT: Sending HL7 results is blocked when an error appears in MeLinDA
- [\[ZH-2183\]](#) - ZH, Daily Work: BlueDiverII - Does not recognize having a pending sample from that machine

Improvement

- [\[ZH-2178\]](#) - Driver TCP Server channel: Limit accepted sockets/connections to 1

New feature

- [\[ZH-2012\]](#) - ZH IFA: Collect and reflect the ZenitPRO warnings regarding some samples

Sprint ZH 2.1.4_Rev9

Bug

- [\[ZH-2147\]](#) - ZH, Configuration, Analyzers, Analyzer / Reagents: A reagent not configured warning is output for an assay not related to the machine
- [\[ZH-2154\]](#) - ZH IFA: Adding a subtest appears without the positivities

Improvement

- [\[ZH-2168\]](#) - Driver HL7, Zenit PRO: Adaptations v1.6.8, change datetime format adding hour minute and second
- [\[ZH-2121\]](#) - ZH, Serumbank, Add tube: Search for the next free position starting from the position occupied by the last added sample
- [\[ZH-2171\]](#) - Driver HL7, ZenitPRO / Fast: Do not process mitosis images
- [\[ZH-2172\]](#) - Driver G-Sight Driver: Receive JP2 images through the WS and not over the network

New feature

- [\[ZH-2170\]](#) - BlueDiverII Driver: Adaptations of the BlueDiver protocol to work with BlueDiverII
- [\[ZH-2025\]](#) - Driver HL7, HOST: Send image analyzer index to LIS